

EPSON RC+ 7.4.1 Release Notes July 11, 2018

Thank you for using EPSON RC+ 7.4.1. This document contains the latest information for this release. Please read before using this software.

DOCUMENTATION	2
GETTING STARTED	2
WHAT'S NEW IN VERSION 7.4.1	2
WHAT'S FIXED IN VERSION 7.4.1	2
VISION GUIDE	3
CAMERA & LENS SELECTION TOOL	3
PARAMETER TUNING TOOL	3



Documentation

All RC+ manuals are in PDF format and are installed on the PC hard disk. These manuals are accessible from the EPSON RC+ 7.0 environment Help menu.

Getting Started

Read the Getting Started chapter in the EPSON RC+ 7.0 User's Guide. This chapter will refer you to the RC700 Controller manual Installation chapter. This contains information for initial connections and start up.

What's New in version 7.4.1

1. Supports N6-850 and N6 clean type robots.

What's Fixed in version 7.4.1

- 1. Added the ability to specify the local coordinate system in Robot Manager Box Wizard.
- 2. Fixed a problem for Robot Manager Pallet Wizard. The arm orientation of the robot was not being recorded when the pallet definition points were taught.
- 3. Fixed a problem for Vision Guide where when Polarity was set to Both, it was displayed as DarkOnLight in the property grid.
- 4. Fixed a problem for Vision Guide where sometimes Correlation could not find a part near the boundary of the search area.
- 5. Fixed a problem for the Timeout property for Geometric and Correlation tools. The tools were not timing out during initial rough search.
- 6. Fixed a problem for Vision Guide Text object UserText property. When the dialog was opened to change the value, the previous value was not being displayed.
- 7. Fixed the VGTuningTool sample. An invalid file was specified for an ImageFile property.
- 8. Fixed IOCondOK result display for Force Guide FollowMove object. The value was not being updated.
- 9. Fixed a problem for Force Guide object execution. When an object with AbortSeqOnFail = True failed during step execution, the objects after the failed object were highlighted.
- 10. Fixed a problem for Force Monitor. If the Remove button is clicked during file remove, an exception occurred.
- 11. Fixed a problem for the Robot Manager Mass Property Wizard. Some labels for the Finish step were incorrect.



Vision Guide

Camera & Lens Selection Tool

A camera and lens selection tool is provided in the EpsonRC70 \Tools folder after installation. A PDF file with instructions is provided in the folder.

Parameter Tuning Tool

A sample project for a vision parameter tuning tool is provided in EpsonRC70\Projects\Samples\Vision\VGTuningTool. A PDF file with instructions is provided in the folder. The GUI Builder option is required to use this tool.